



Creating Greater Certainty in the
Fight Against Cancer

Poster No. 3003
AACR 2021


In Vitro and In Vivo Model for Evaluation of Novel Therapeutic Agents in Patient derived- Lymphoblastic Leukemia (B-ALL)

Bandana Ajay Vishwakarma, PhD, Bhavna Verma, PhD, and
Amy Wesa, PhD

Champions Oncology, Rockville, MD

B Cell Acute Lymphoblastic Leukemia (B-ALL)

- Malignant transformation and proliferation of lymphoid progenitors in peripheral blood, bone marrow, and extramedullary sites.
- Prevalent in children and young adults
- 6150 new cases and 1520 deaths
- Relapse and resistant cases
- Drug screening platform - ex vivo and in vivo

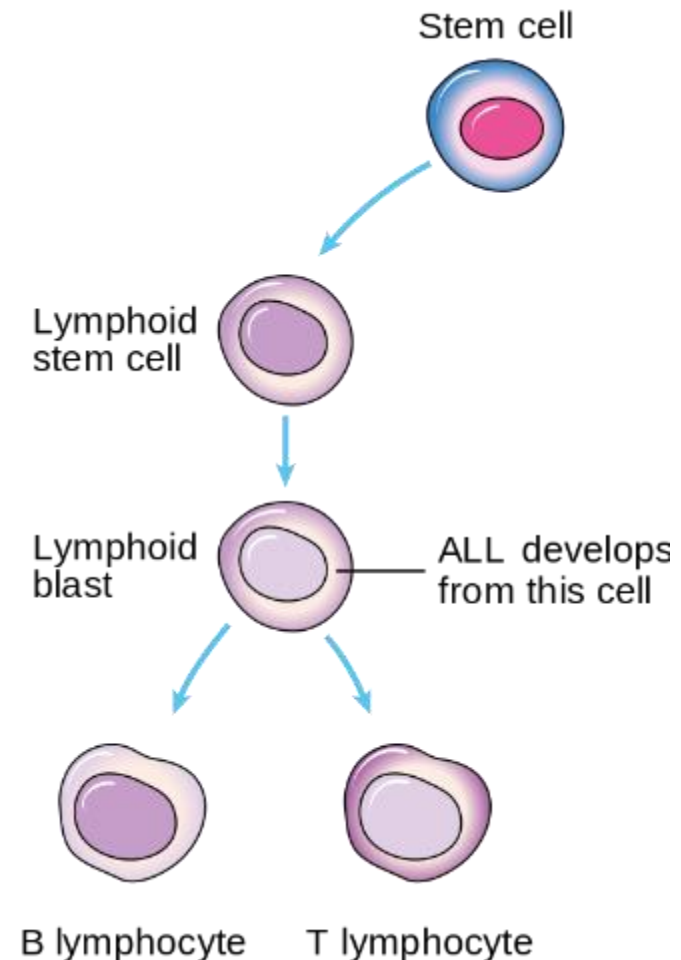

Champions B-ALL Model Characterization


- **Clinical Characteristics:**
 - Age, Gender, Race
- **Precursor B Cell ALL**
- **De Novo**
- **MYC, ATM, NRAS, PTPN11, RUNX1**
- **Immunophenotyped**
- **t(9;22) [BCR-ABL]+: 13 models**
- **t(9;22) [BCR-ABL]-: 11 models**

Champions Models	Sample Type	Status Collection	Cytogenetics Abnormality	Clinical Flow Cytometry Phenotype
CTG-2406	Pheresis	Unknown	N/A	N/A
CTG-2447	Pheresis	Unknown	N/A	CD10 bright +CD19+CD20 dim, variable, small subset + CD34 dim, variable, small subset +CD79+HLA-DR bright +TdT+
CTG-2409	Pheresis	De Novo	42,XX,+8,t(9;22)(q34;q11.2),+der(22)t(9;22)[1]/46,XX[1]	CD5-CD10+CD13-CD14-CD15-CD19+CD20 var +CD33-CD34 var +CD56-CD64-CD79 a +HLA-DR+TdT var +
CTG-3682	PB-MNC	De Novo	42,XX,+8,t(9;22)(q34;q11.2),+der(22)t(9;22)[1]/46,XX[1]	CD5-CD10+CD13-CD14-CD15-CD19+CD20 var +CD33-CD34 var +CD56-CD64-CD79 a +HLA-DR+TdT var +
CTG-3683	PB-MNC	De Novo	46,XY[5] (Normal)	CD5-CD10+CD13-CD14-CD15-CD19+CD20dim/variable +CD33CD34+CD56CD64CD79+HLADR+TdTdim/variable +
CTG-3684	PB-MNC	De Novo	Normal	CD5-CD10 var +CD13 var +CD14-CD15-CD19 var+ CD20-CD33 var+CD34+CD38 var +CD56-CD64-CD79 var +HLA-DR+TdT+
CTG-2707	Pheresis	De Novo	45,XY,-7,t(9;22)(q34;q11.2)[20]	CD5-CD10+CD13+CD14-CD15-CD19+CD20-CD33 var + CD34 var +CD56-CD64-CD79+HLA-DR bright +TdT+
CTG-3685	Pheresis	De Novo	46,XX,t(4;11)(q21;q23)[6]/46,XX[5]	CD5-CD10-CD13-CD14-CD15 aberrant expression + CD19+CD20-CD33-CD34-CD38+CD56-CD64-CD79+ HLA-DR+TdT+
CTG-3686	PB-MNC	De Novo	47,XY,del(7)(p15),+add(8)(p11.2),der(9)del(9)(p13)t(9;22)(q34;q11.2),der(22)t(9;22)[7]/ 46,XY[1]	CD5-CD10+CD13-CD14-CD15-CD19+CD20 var +CD33-CD34+CD38-CD56-CD64-CD79+HLA-DR+TdT+

Development of ex vivo and in vivo platform for evaluation of novel therapeutics in patient-derived B-ALL

Schematic Representation of the Ex Vivo Study

Patient-derived B-ALL models

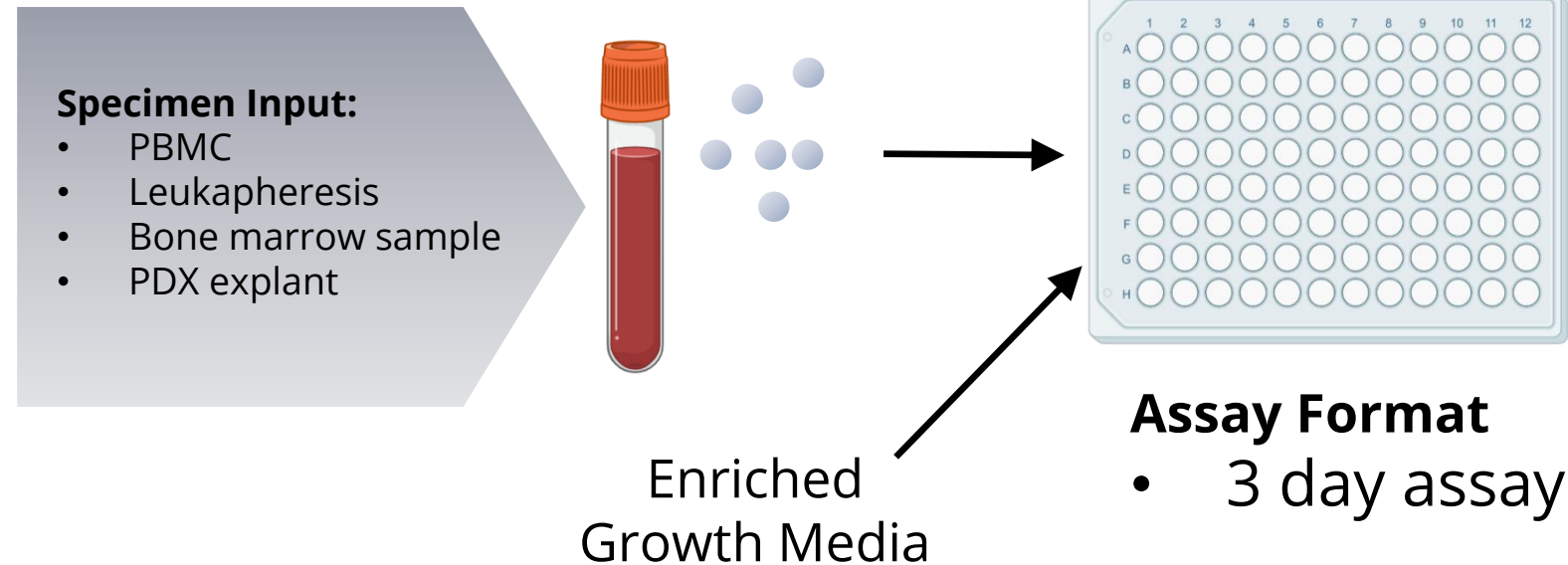


Applications

- Drug Screen
- Biomarker Discovery
- Mechanism of Action
- In vivo Model Selection

Study Endpoints:


- Proliferation
- Tumor Cytotoxicity
- Apoptosis
- Phenotyping

Evaluation of Culture Conditions for 3 day B-ALL cultures

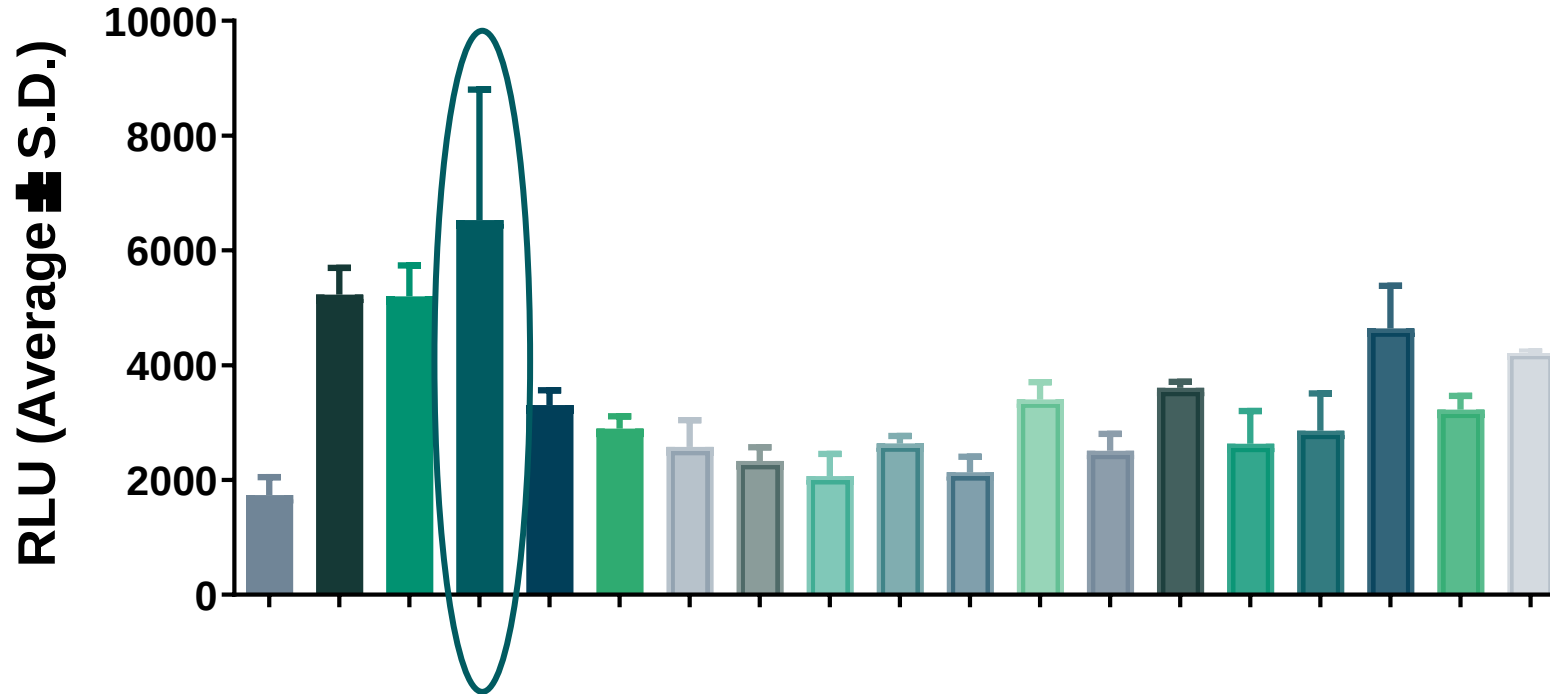


Selection of media, supplements and other culture conditions (representative of several studies with similar results across models)





Reproducible Performance in Proliferation/ Survival in Enriched Media

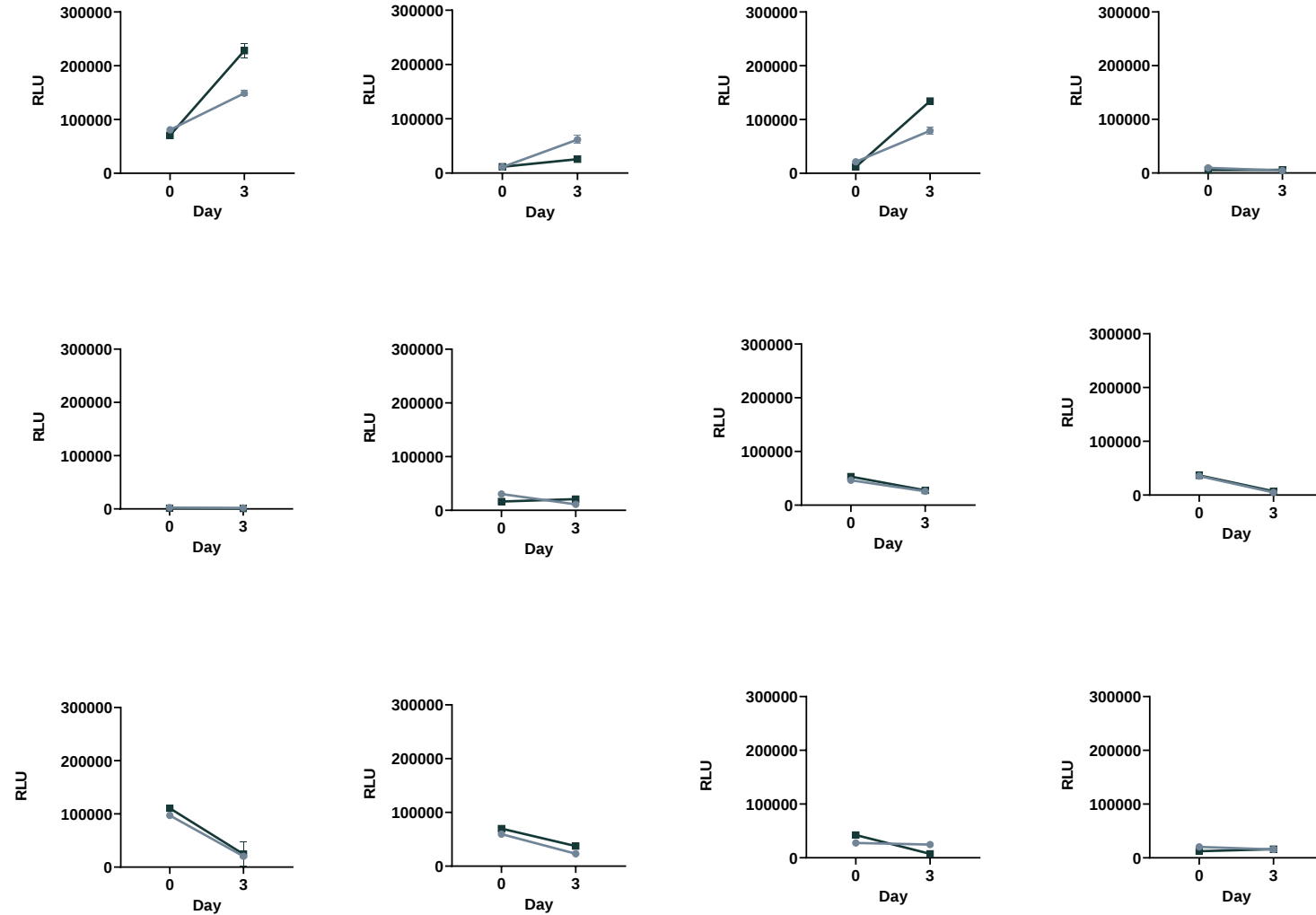



Models Maintain B-ALL Phenotype in Culture

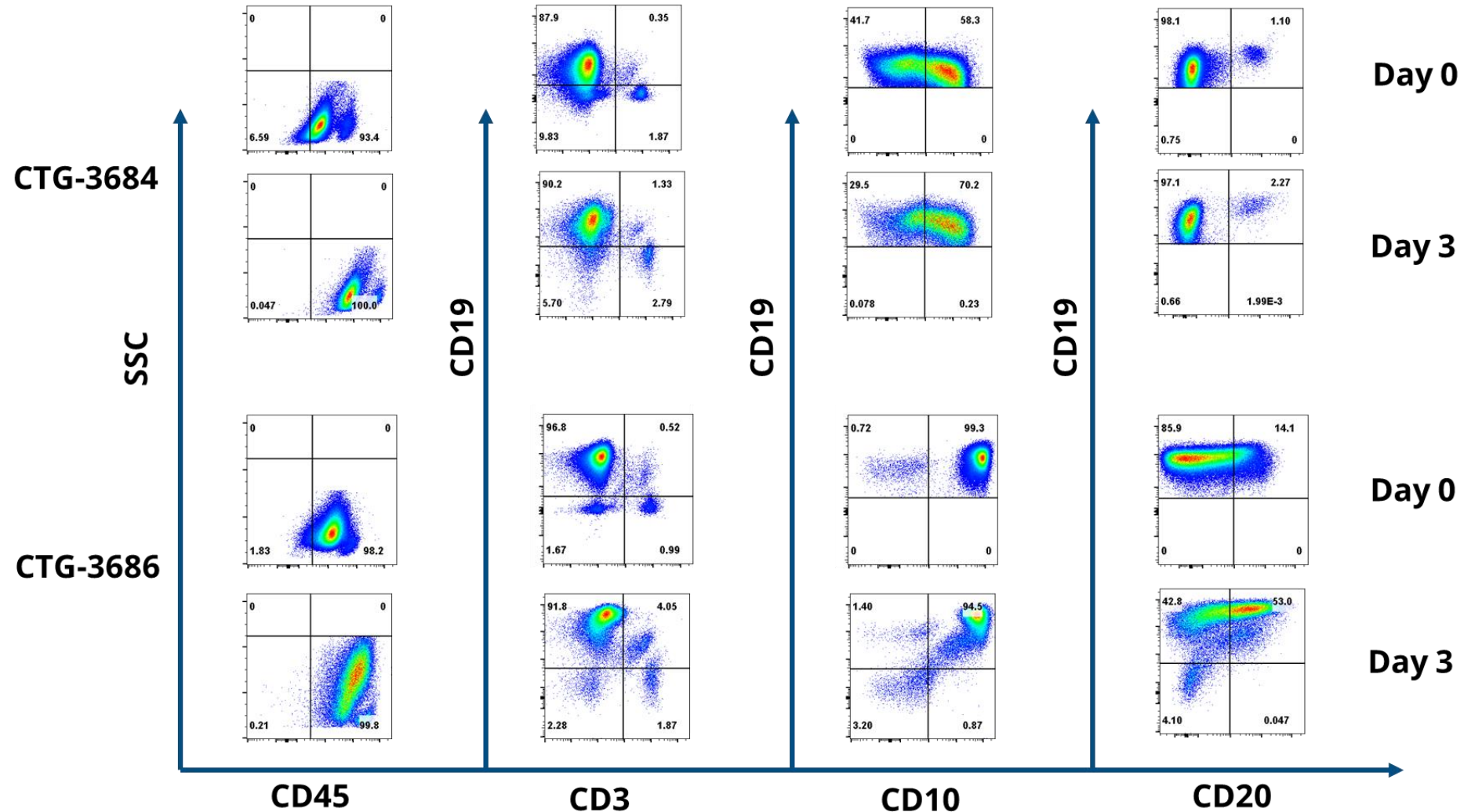

Responsiveness of B-ALL to Standard of Care Treatment in Ex Vivo Platform

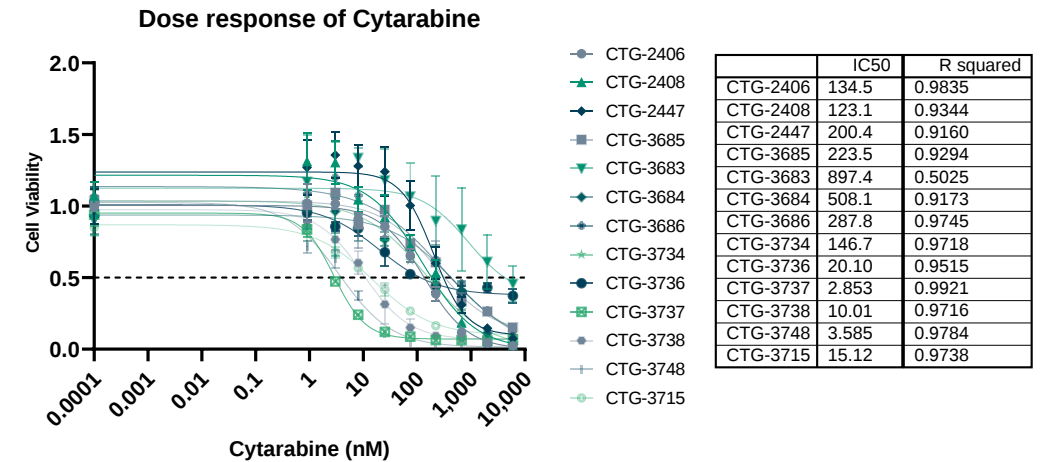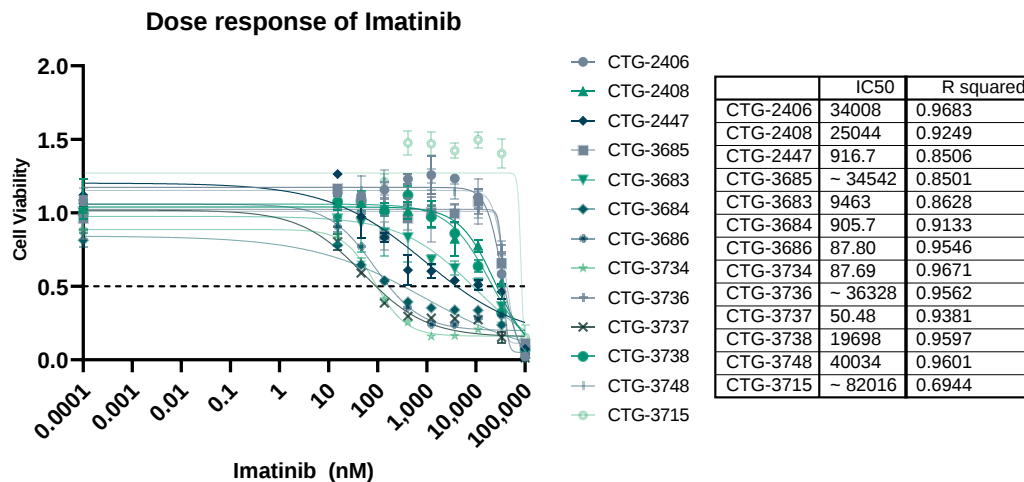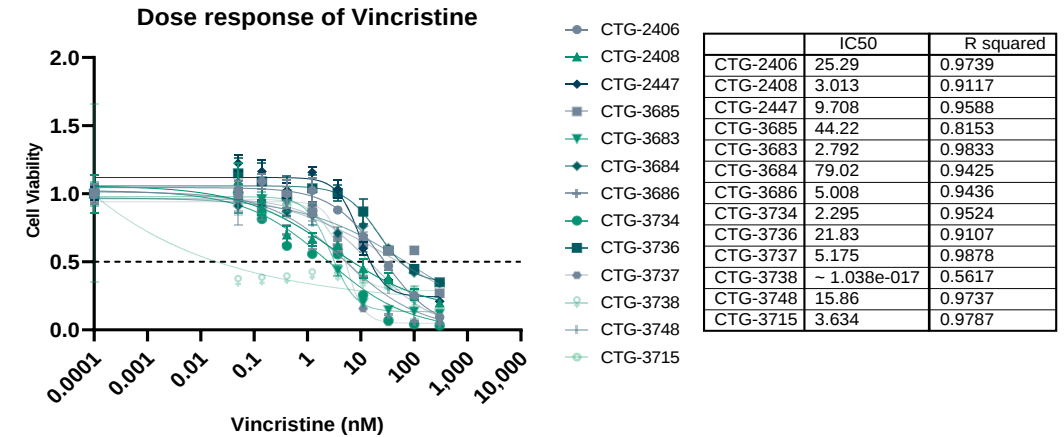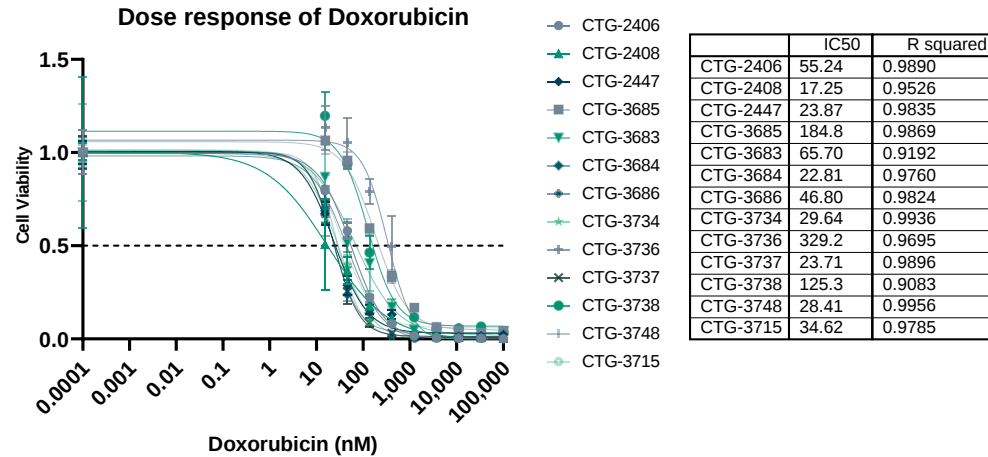


Mutational Status Drives Differential Response to Imatinib


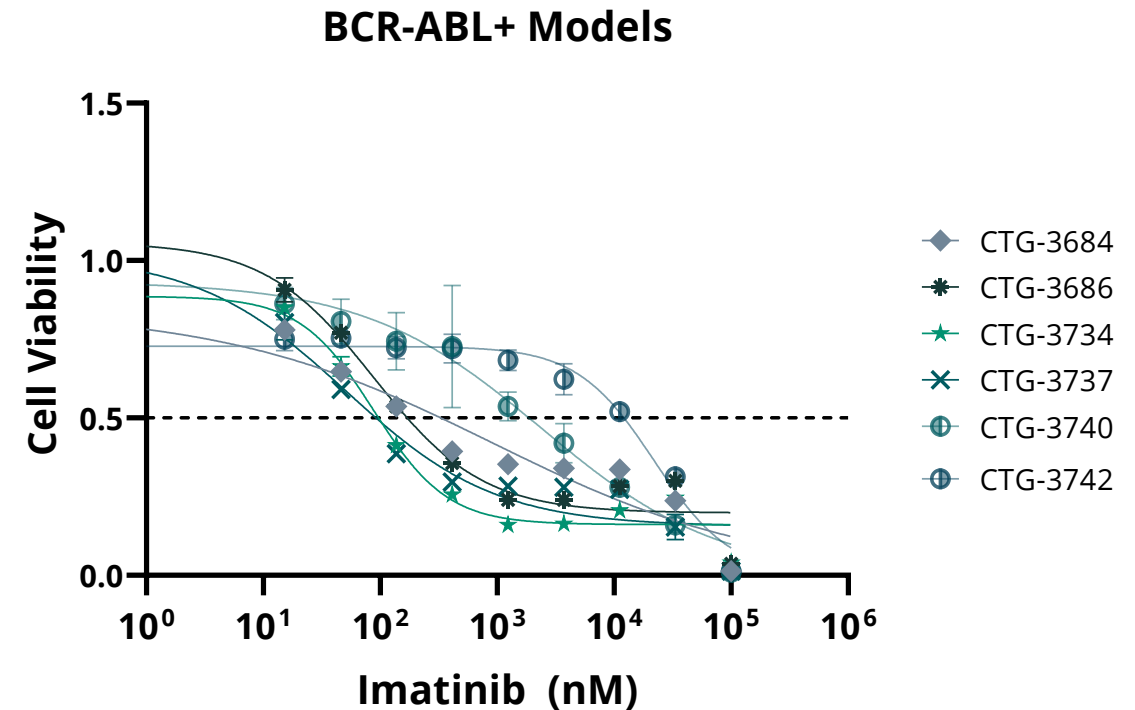
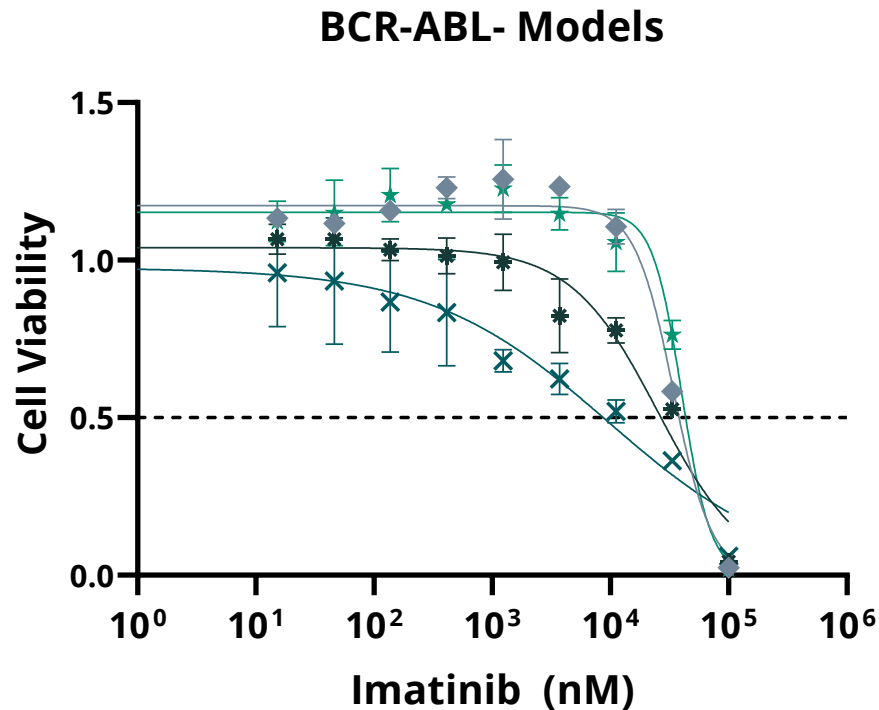


Champions' B-ALL PDX Development for In Vivo Drug Efficacy Study

Engraftment and Kinetics of Primary B-ALL cells In Vivo


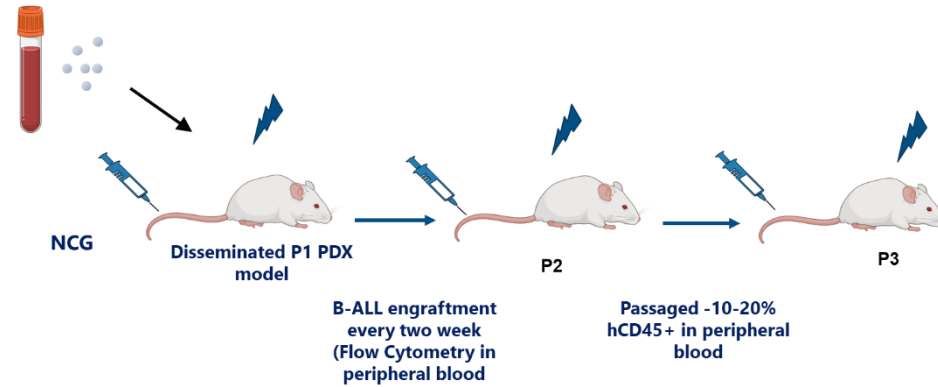


Schematic representation of B-ALL PDX development

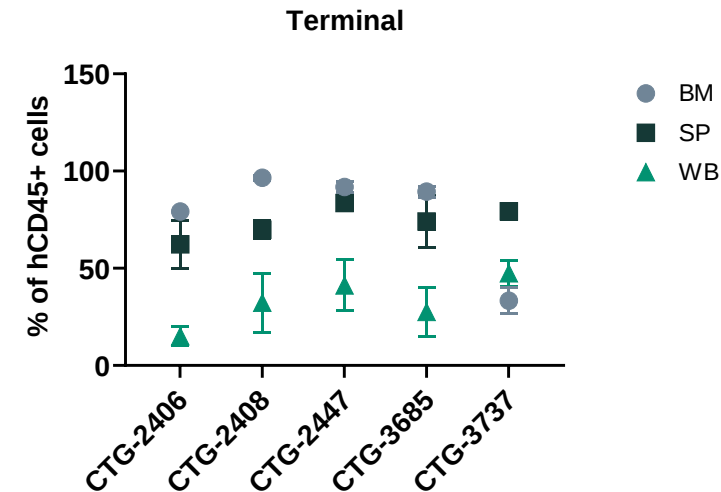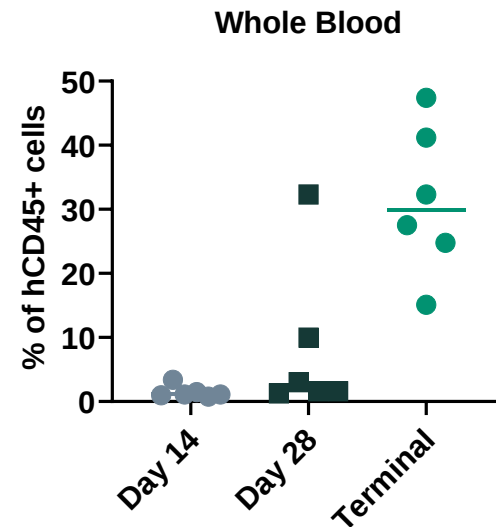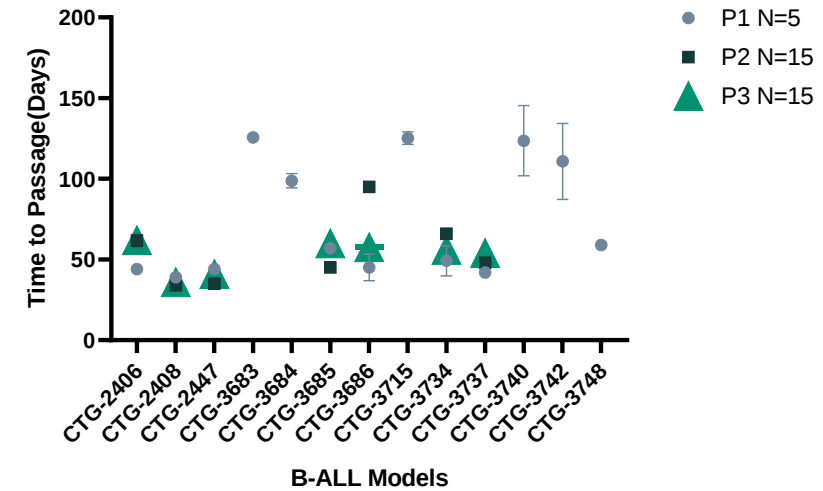

Immunophenotype of Patient Sample is Comparable to P1 BM and Spleen Cells




CTG-2408

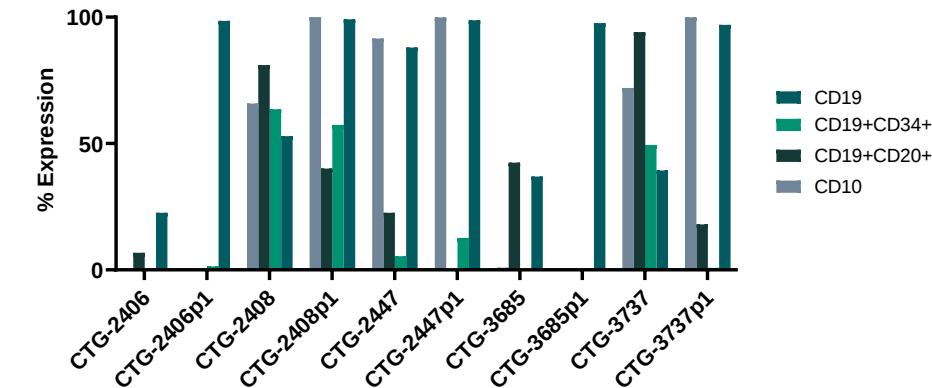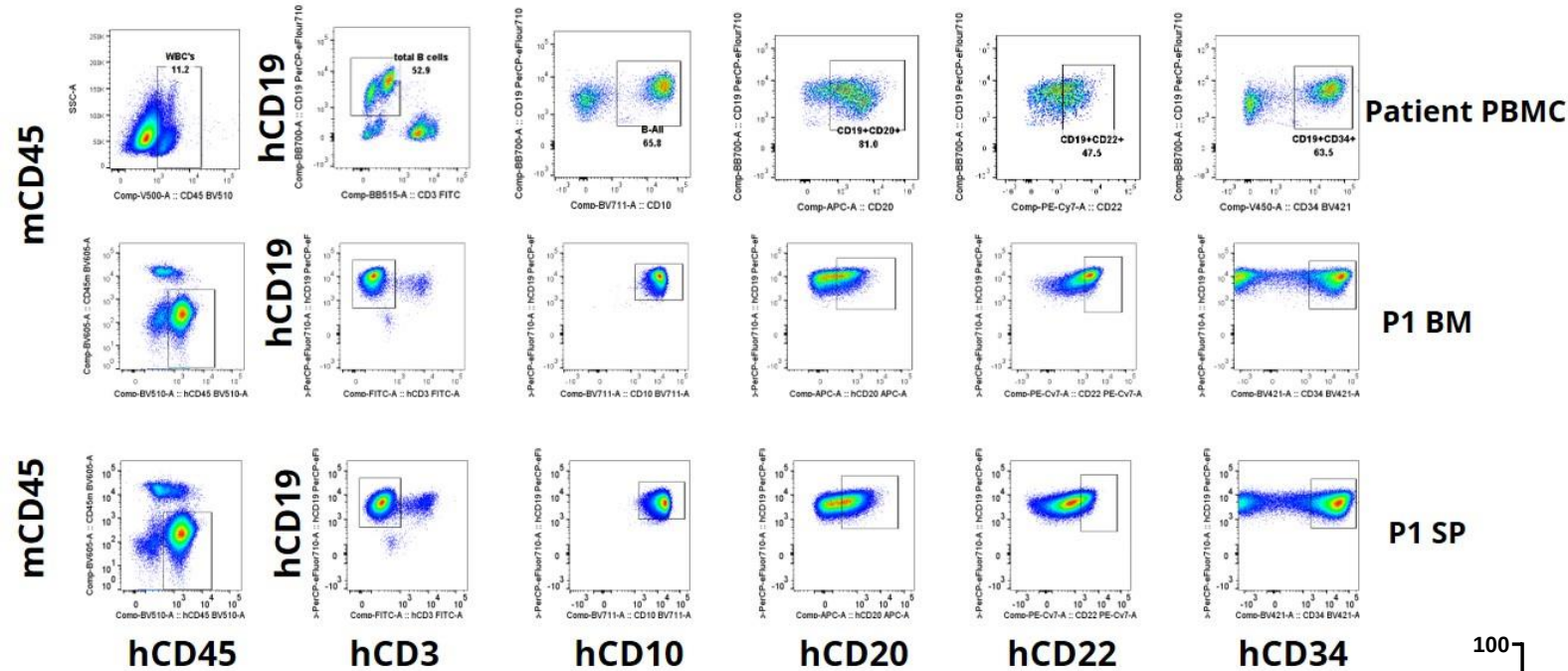

Comparable Ex Vivo Response of Primary B-ALL and Passaged B-ALL PDX (P2) Models







	Model A CTG-2406		Model B CTG-2408		Model C CTG-2447	
SOC	Patient Specimen	B-ALL PDX Explant (Splenocytes)	Patient Specimen	B-ALL PDX Explant (Splenocytes)	Patient Specimen	B-ALL PDX Explant (Splenocytes)
Doxorubicin	IC50 – 32.28nM R ² -0.99	IC50 – 23.82nM R ² -0.99	IC50 – 17.25nM R ² -0.95	IC50 – 0.56nM R ² -0.95	IC50 – 23.87nM R ² -0.98	IC50 – 15.09nM R ² -0.99
Vincristine	IC50 – 20.27nM R ² -0.96	IC50 – 6.97nM R ² -0.96	IC50 – 3.013nM R ² -0.91	IC50 – 1.579nM R ² -0.84	IC50 – 9.708nM R ² -0.95	IC50 – 31.77nM R ² -0.98
Cytarabine	IC50 – 828.4nM R ² -0.95	IC50 – 828.8nM R ² -0.90	IC50 – 123.1nM R ² -0.93	IC50 – 473.7nM R ² -0.88	IC50 – 200.4nM R ² -0.916	IC50 – 83.85nM R ² -0.96
Imatinib	IC50 – 21.16uM R ² -0.94	IC50 – 30.10uM R ² -0.93	IC50 – 25.04uM R ² -0.92	No Response	IC50 – 0.916uM R ² -0.85	IC50 – 3.755uM R ² -0.85

Venetoclax Treatment Inhibits Proliferation of Patient-derived B-ALL Cells (CTG-2408) In Vivo




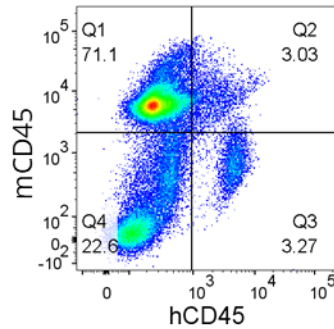
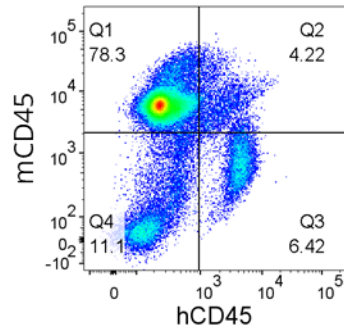
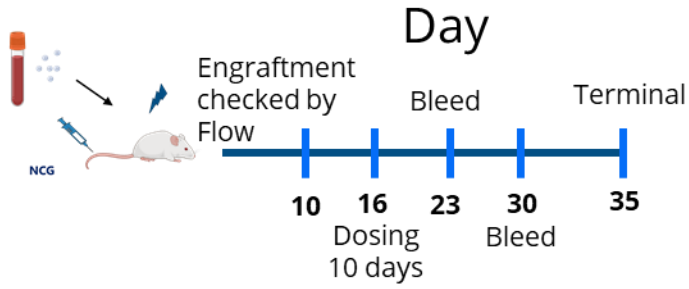


Bone Marrow Day 10 before dosing

Vehicle Treated

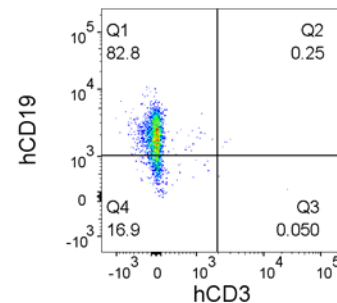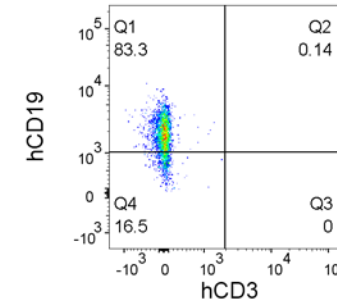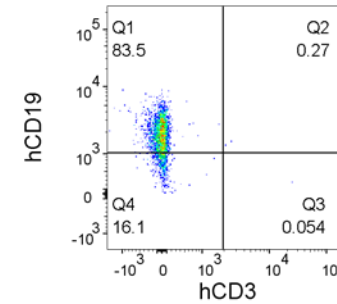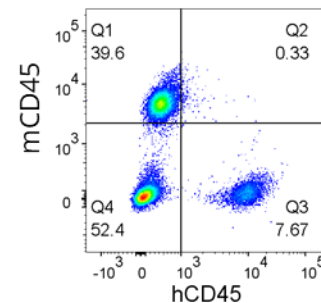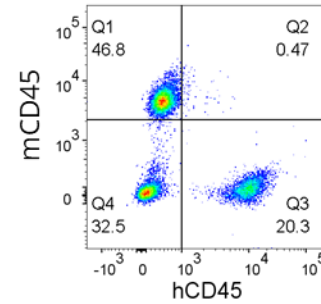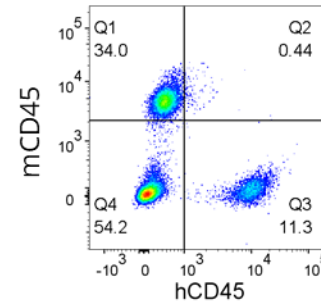

Venetoclax Treated

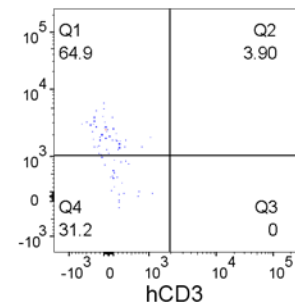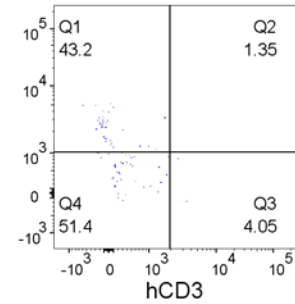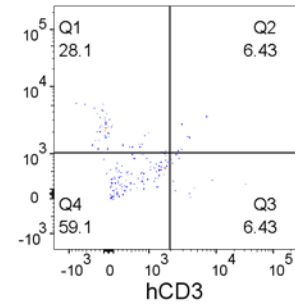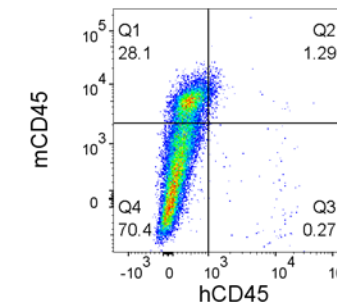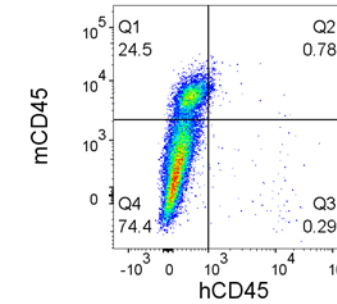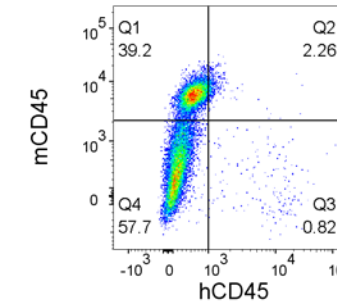

Peripheral Blood Day 7 of treatment

Summary & Next Steps

- Patient derived primary B-ALL cells survive and proliferate in enriched media.
- B-ALL models showed responsiveness to standard of care drugs in ex vivo platform. Interestingly, molecular subtyping (BCR-ABL+) demonstrated correspondence with drug response to imatinib in this platform.
- NCG mice support the engraftment and development of B-ALL disease initiating and leukemic cells in vivo.
- The passaged patient-derived preclinical xenograft model recapitulates the B-ALL disease.
- Venetoclax inhibits the proliferation of patient derived B-ALL cells in vivo suggesting that the patient-derived preclinical in vivo model can be used for drug efficacy studies.
- Passaged B-ALL models will provide a renewable source for both ex vivo and in vivo platforms for drug screening/drug discovery.
- WES and RNA-Seq is in progress for all models.



Creating Greater Certainty in the
Fight Against Cancer



www.ChampionsOncology.com

Corporate Headquarters

One University Plaza, Suite 307
Hackensack, NJ 07601

Research US Headquarters

1330 Piccard Drive, Suite 025
Rockville, MD 20850

Research EU Facilities

260, Avenue Marie Curie
74160 Archamps, France